



OOSAFE® IVF LABORATORY DISINFECTANTS

used by leading IVF clinics in more than 100 countries

Manufactured in
an EN ISO 13485:2016
certified facility.



GOosafe in IVF Laboratories

OOSAFE® DISINFECTANT FOR CO₂ INCUBATORS AND LAMINAR FLOW HOODS FOR IVF LABORATORY



High quality

- Ready-to-use solution
- Anticorrosive
- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection/low evaporation rate
- Fast disinfection effect

Health and safety

- MEA and HSSA tested
- Transparent
- Odorless
- Alcohol-free
- No VOC release


Excellent material compatibility

- Glass, plastic, acrylic and metal

Directions:

Spray: Apply undiluted, spray and distribute evenly on the surface with a sterile cloth. Wait 15 minutes for the disinfection effect.

Wipes: Wipe the surface evenly and wait 15 minutes for the disinfection to take effect.

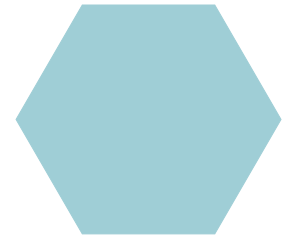

BACTERICIDAL

- 15 min Pseudomonas aeruginosa
15 min Staphylococcus aureus
15 min Enterococcus hirae

FUNGICIDAL

- 15 min Candida albicans

MYCOBACTERICIDAL AND TUBERCULOCIDAL

- 15 min Mycobacterium terrae
15 min Mycobacterium avium

VIROCIDAL

- 15 min Vaccinia virus*
15 min Poliovirus
5 min Adenovirus

*Effective against all enveloped viruses including HIV, Hepatitis B and C, Herpes simplex, Influenza A, Ebola

TIME UNTIL FULL EFFECT

MICROORGANISM



Product code:
OODW-70
Oosafe® Disinfection - Wipes,
70 wipes per bucket.



Product code:
OODIH-1000
Oosafe® Disinfectant for CO₂
Incubators and Laminar Flow Hoods,
1 Liter bottle with spray



Product code:
OODIH-5000
Oosafe® Disinfectant for CO₂
Incubators and Laminar Flow Hoods,
5 Liters refill container



Product code:
OODIH-0050
Oosafe® Disinfectant for CO₂
Incubators and Laminar Flow Hoods,
50 ml in spray bottle

OOSAFE® HAND DISINFECTANT FOR IVF LABORATORY



Health and safety

- MEA and HSSA tested
- Transparent
- Odorless
- Alcohol-free
- No VOC release
- Keeps your skin soft

High quality

- Ready-to-use solution
- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection/low evaporation rate
- Fast disinfection effect

Directions:

Follow laboratory's hygienic procedure for washing hands before using hand disinfectant. For effective hand disinfection rub hands, apply minimum 2 pumps (3 ml) of disinfectant and rub hands till they are completely dry.

TIME UNTIL FULL EFFECT

BACTERICIDAL

Pseudomonas aeruginosa

30 sec
Staphylococcus aureus
Enterococcus hirae

Escherichia coli

FUNGICIDAL

30 sec
Candida albicans

VIRUCIDAL

30 sec
Vaccinia virus*

MICROORGANISM

**Effective against all enveloped viruses including HIV, Hepatitis B and C, Herpes simplex, Influenza A, Ebola*



Product code:
OODH-0500
Oosafe® Hand Disinfectant,
500 ml bottle with pump



Product code:
OODH-5000
Oosafe® Hand Disinfectant,
5 Liters refill container



Product code:
OODH-0050
Oosafe® Hand Disinfectant,
50 ml in spray bottle

OOSAFE® LABORATORY SURFACE DISINFECTANT FOR IVF LABORATORY



Health and safety

- MEA and HSSA tested
- Clear - slightly yellow
- Alcohol-free
- Very low level of VOC release

High quality

- Securely sealed
- Valid for 3 years, even after opening
- Long term disinfection

Excellent material compatibility

- Laboratory surfaces and floors
- Glass, Plastic, Tiles, Rubber

Directions: Dilute with water (according to the table below) and apply diluted cleaning agent on laboratory floors and surfaces. After application wait 30 minutes for the required disinfection effect and rinse the surface with water.

MICROORGANISM

BACTERICIDAL

TIME UNTIL FULL EFFECT	MICROORGANISM	DILUTION
15 min	Pseudomonas aeruginosa	1:50
	Staphylococcus aureus	1:50
	Enterococcus hirae	1:50

FUNGICIDAL

15 min	Candida albicans	1:50
--------	------------------	------

MYCOBACTERICIDAL AND TUBERCULOCIDAL

30 min	Mycobacterium terrae	1:25
	Mycobacterium avium	1:25

VIRUCIDAL

15 min	Vaccinia virus*	1:50
30 min	Poliovirus type 1	1:25
10 min	Adenovirus type 5	1:25

*Effective against all enveloped viruses including HIV, Hepatitis B and C, Herpes simplex, Influenza A, Ebola



Product code:
OOSDF-02000
Oosafe® Laboratory Surface Disinfectant, 2 Liters bottle with dose cup



Product code:
OOSDF-10000
Oosafe® Laboratory Surface Disinfectant, 10 Liters refill container



Product code:
OOSDF-00050
Oosafe® Laboratory Surface Disinfectant, 50 ml in bottle



we support



THE DANISH CHILDHOOD
CANCER ORGANISATION



 facebook.com/sparMED

 twitter.com/sparMED

 linkedin.com/company/sparmed-aps